


APPLICATIONS

Extraction and Analysis of Fentanyl and Analogs from Whole Blood Using a Kinetex[®] F5 LC Column

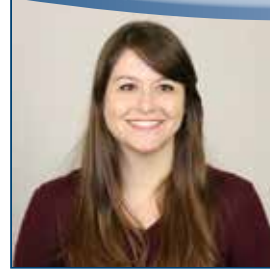

Charissa Goggin
Technical Sales Consultant

In addition to being a chromatography enthusiast, Charissa enjoys home improvement projects and has renovated an entire house with her family!

Charissa Goggin¹ and Szabolcs Sofalvi²

¹Phenomenex Inc., 411 Madrid Ave, Torrance, CA 90501 USA

²Cuyahoga County Medical Examiner's Office, Cuyahoga County Regional Forensic Science Laboratory, 11001 Cedar Avenue, Cleveland, OH 44106 USA

Introduction

Fentanyl and fentanyl analogs are high potency synthetic opioids that are 80-100x more powerful than morphine. Although fentanyl use is strictly regulated, new fentanyl analogs keep emerging with molecular structures that are closely related, making separation, identification, and quantitation difficult for analysts. This method explores the identification of 26 compounds and quantitation of 24 compounds using multiple sample preparation techniques to clean-up the whole blood matrix. A core-shell LC column was utilized for its high efficiency and selective properties and the F5 phase was chosen for separating geometric and structural isomers.

Methods

Sample Pre-treatment

500 µL of a whole blood sample was lysed osmotically with 500 µL of DI Water and internal standard was added. The samples were vortexed for 7 seconds and they sat for 10 minutes. Protein precipitation was accomplished by drop-wise addition of 1.5 mL of ice cold methanol/acetonitrile (10:90) while vortexing. The samples were then incubated for 10 min in the freezer and then centrifuged at 3,500 rpm (2,800 xg) for 10 min. The supernatants were poured into new 16 x 125 mm test tubes followed by the addition of 3 mL DI water and vortexed for 7 sec.

SPE Conditions

- Cartridge:** Strata[®] Screen C (55 µm, 70 Å), 200 mg/6 mL
- Part No.:** 8B-S016-FCH
- Condition:** 3 mL of Methanol
- Equilibrate:** 2 mL of DI water
- Load:** Pre-treated samples were loaded onto the cartridge under 2 psi
- Wash 1:** 2 mL of DI water (All 3 wash steps were eluted under 5 psi positive pressure)
- Wash 2:** 1.5 mL of 1.0 M Acetic acid
- Wash 3:** 2 mL of Methanol
- Dry:** Cartridge for 10 minutes
- Elute:** 2x 1.5 mL of Dichloromethane/2-propanol/Ammonium hydroxide (78:20:2) under 5 psi. Full flow was applied at the end of each elution step for 30 sec.
- Dry Down:** Evaporated to dryness
- Reconstitute:** 40 µL of initial mobile phase. The samples were then vortexed for 7 sec and centrifuged at 2,800 xg for 5 min. The supernatants were transferred to conical inserts for UHPLC-MS/MS analysis.

LC Conditions


- Column:** Kinetex[®] 2.6 µm F5
- Dimensions:** 150 x 2.1 mm
- Part No.:** 00F-4723-AN
- Mobile Phase:** A: 0.1% Formic acid in Water
B: 0.1% Formic acid in Acetonitrile
- Gradient:**

Time (min)	% B
0	20
17	25
17.25	95
18.75	95
19.5	20
21	20
- Flow Rate:** 0.5 mL/min
- Column Oven Temperature:** 40 °C
- Autosampler Temperature:** 5 °C
- Detector:** MS/MS (Thermo TSQ Quantis[™] Triple Quad Tandem)

	Analyte Name	Retention Time (min)
1.	Norfentanyl	1.19
2.	N-methyl norfentanyl	1.19
3.	Norcarfentanil	1.47
4.	Methoxyacetyl fentanyl	3.07
5.	Acetyl fentanyl	3.51
6.	β-hydroxy fentanyl	3.67
7.	Benzyl fentanyl	3.95
8.	para-methoxyacetyl fentanyl	4.67
9.	Alfentanil	5.02
10.	4-ANPP	5.50
11.	Acryl fentanyl	5.93
12.	Fentanyl	6.42
13.	para-fluoroacryl fentanyl	8.01
14.	Cyclopropyl fentanyl	8.59
15.	para-fluorofentanyl	8.82
16.	2-furanyl fentanyl	9.04
17.	(E)-crotonyl fentanyl	9.35
18.	(±)-trans-3-methylfentanyl	9.74
19.	Fentanyl carbamate	10.27
20.	(±)-cis-3-methylfentanyl	10.31
21.	Carfentanil	11.40
22.	Butyryl fentanyl	11.66
23.	Sufentanil	13.76
24.	para-fluoroisobutyryl fentanyl	14.29
25.	Phenyl fentanyl	15.13
26.	Cyclopentenyl fentanyl	15.59

Results

Figure 1.
Separation and Analysis of Fentanyl and Metabolites Using Kinetex[®] F5 LC Column

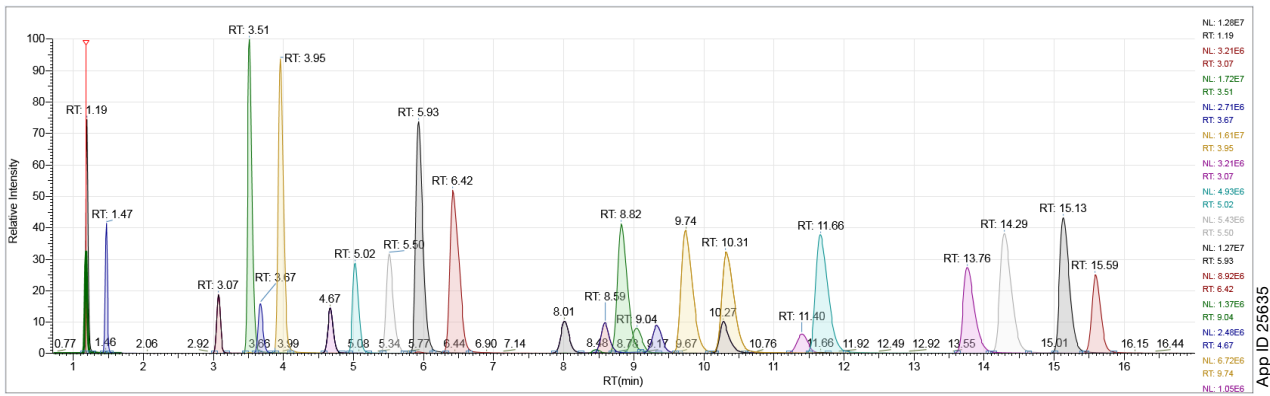

Conclusion

Using whole blood as a matrix requires pre-treatment and sample preparation to clean-up prior to analysis. Employing a solid phase extraction method allows for matrix clean-up and the sample is concentrated which leads to more sensitive LC-MS/MS results. The Kinetex F5 LC column utilized a core-shell media to help which has 5 modes of selectivity: hydrophobic, aromatic, electrostatic, steric, and hydrogen bonding. This makes the phase extremely useful for isomeric analysis, and a great choice for fentanyl analog analysis.

References

1. Szabolcs Sofalvi, Eric S Lavins, Ian T Brooker, Claire K Kaspar, John Kucmanic, Carrie D Mazzola, Christie L Mitchell-Mata, Cassandra L Clyde, Rindi N Rico, Luigino G Apollonio, Charissa Goggin, Brittany Marshall, Danielle Moore, Thomas P Gilson, (2019). Unique Structural/Stereo-Isomer and Isobar Analysis of Novel Fentanyl Analogues in Postmortem and DUID Whole Blood by UHPLC-MS-MS, *Journal of Analytical Toxicology*, Volume 43, Issue 9, bkz056, <https://doi.org/10.1093/jat/bkz056>

Phenomenex would like to thank Szabi Sofalvi and the rest of the Toxicology Department at Cuyahoga County Medical Examiner's Office for the use of their data and their time in developing this method.

Ordering Information

Kinetex® LC Columns

5 µm Minibore Columns (mm)				SecurityGuard™ ULTRA Cartridges†
Phases	50 x 2.1	100 x 2.1	150 x 2.1	3/pk
F5	00B-4724-AN	00D-4724-AN	00F-4724-AN	AJO-9322 for 2.1 mm ID

5 µm MidBore™ Columns (mm)			SecurityGuard ULTRA Cartridges†
Phases	100 x 3.0	150 x 3.0	3/pk
F5	00D-4724-YO	00F-4724-YO	AJO-9321 for 3.0 mm ID

5 µm Analytical Columns (mm)					SecurityGuard ULTRA Cartridges†
Phases	50 x 4.6	100 x 4.6	150 x 4.6	250 x 4.6	3/pk
F5	00B-4724-E0	00D-4724-E0	00F-4724-E0	00G-4724-E0	AJO-9320 for 4.6 mm ID

2.6 µm Minibore Columns (mm)					SecurityGuard ULTRA Cartridges†
Phases	30 x 2.1	50 x 2.1	100 x 2.1	150 x 2.1	3/pk
F5	00A-4723-AN	00B-4723-AN	00D-4723-AN	00F-4723-AN	AJO-9322 for 2.1 mm ID


2.6 µm MidBore™ Columns (mm)				SecurityGuard ULTRA Cartridges†
Phases	50 x 3.0	100 x 3.0	150 x 3.0	3/pk
F5	00B-4723-Y0	00D-4723-Y0	00F-4723-Y0	AJO-9321 for 3.0 mm ID


2.6 µm Analytical Columns (mm)				SecurityGuard ULTRA Cartridges†
Phases	50 x 4.6	100 x 4.6	150 x 4.6	3/pk
F5	00B-4723-E0	00D-4723-E0	00F-4723-E0	AJO-9320 for 4.6 mm ID

1.7 µm Minibore Columns (mm)				SecurityGuard ULTRA Cartridges†
Phases	50 x 2.1	100 x 2.1	150 x 2.1	3/pk
F5	00B-4722-AN	00D-4722-AN	00F-4722-AN	AJO-9322 for 2.1 mm ID

† SecurityGuard ULTRA Cartridges require holder, Part No.: AJO-9000.

Strata® Screen C SPE

Format	Sorbent Mass	Part Number	Unit
Tube			
	100 mg	8B-S016-EAK	1 mL (100/box)
	100 mg	8B-S016-EBJ	3 mL (50/box)
	150 mg	8B-S016-SBJ	3 mL (50/box)
	150 mg	8B-S016-SCH	6 mL (30/box)
	200 mg	8B-S016-FBJ	3 mL (50/box)
	200 mg	8B-S016-FCH	6 mL (30/box)
	300 mg	8B-S016-RBJ	3 mL (50/box)
	300 mg	8B-S016-RCH	6 mL (30/box)
	500 mg	8B-S016-HCH	6 mL (30/box)

96-Well Plate			
	50 mg	8E-S016-DGB	2 Plates/Box

APPLICATIONS

Australia

t: +61 (0)2-9428-6444
auinfo@phenomenex.com

Austria

t: +43 (0)1-319-1301
anfrage@phenomenex.com

Belgium

t: +32 (0)2 503 4015 (French)
t: +32 (0)2 511 8666 (Dutch)
beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681
info@phenomenex.com

China

t: +86 400-606-8099
cninfo@phenomenex.com

Denmark

t: +45 4824 8048
nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063
nordicinfo@phenomenex.com

France

t: +33 (0)1 30 09 21 10
franceinfo@phenomenex.com

Germany

t: +49 (0)6021-58830-0
anfrage@phenomenex.com

India

t: +91 (0)40-3012 2400
indiainfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405
eireinfo@phenomenex.com

Italy

t: +39 051 6327511
italiainfo@phenomenex.com

Luxembourg

t: +31 (0)30-2418700
nlinfo@phenomenex.com

Mexico

t: 01-800-844-5226
tecnicomx@phenomenex.com

The Netherlands

t: +31 (0)30-2418700
nlinfo@phenomenex.com

New Zealand

t: +64 (0)9-4780951
nzinfo@phenomenex.com

Norway

t: +47 810 02 005
nordicinfo@phenomenex.com

Portugal

t: +351 221 450 488
ptinfo@phenomenex.com

Singapore

t: +65 800-852-3944
sginfo@phenomenex.com

Spain

t: +34 91-413-8613
esinfo@phenomenex.com

Sweden

t: +46 (0)8 611 6950
nordicinfo@phenomenex.com

Switzerland

t: +41 (0)61 692 20 20
swissinfo@phenomenex.com

Taiwan

t: +886 (0) 0801-49-1246
twinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367
ukinfo@phenomenex.com

USA

t: +1 (310) 212-0555
info@phenomenex.com

**☉ All other countries/regions
Corporate Office USA**

t: +1 (310) 212-0555
info@phenomenex.com

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at international@phenomenex.com



Your happiness is our mission. Take 45 days to try our products. If you are not happy, we'll make it right.

www.phenomenex.com/behappy

Terms and Conditions

Subject to Phenomenex Standard Terms and Conditions which may be viewed at www.phenomenex.com/TermsAndConditions.

Trademarks

Strata and Kinetex are registered trademarks and SecurityGuard, Be-Happy, and Mid-Bore are trademarks of Phenomenex. TSQ Quantis is a trademark of Thermo Fisher Scientific Inc.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures.

© 2019 Phenomenex, Inc. All rights reserved.