


Circumference	Cap Type
11.8 cm	GL38
12.6 cm	GL40/S40
14.1 cm	GL45
16.8 cm	DIN51
18.5 cm	DIN61

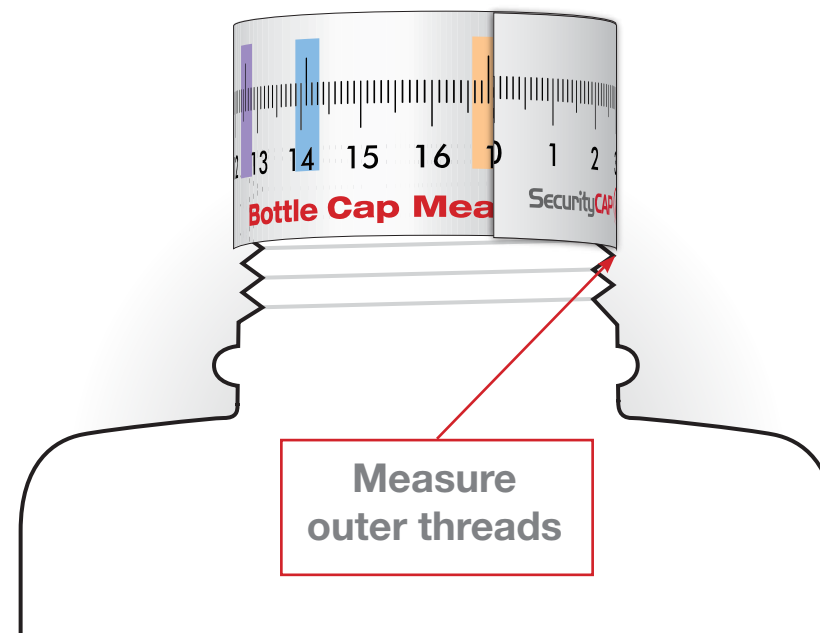

Print

1. Download and print.pdf
2. Be sure to select “actual size” or scale = 100%
3. Cut along the dotted line

Use

1. Remove bottle cap.
2. Place the end of the measuring tape (20 cm side) on your solvent bottle or waster container at the threads.
3. Wrap the tape around the threads until the beginning (0 cm side) aligns with the rest of tape measure.
4. Read indicating color to determine correct Cap Type*

* If tape does not line up with a color indicator please contact your Phenomenex representative.