



## TN-1321

# Next-level Sensitivity for the Quantification of Warfarin and Furosemide in Human Plasma

Ian Moore<sup>1</sup>, Rahul Baghla<sup>2</sup>, and Bryan Tackett, PhD<sup>3</sup>

<sup>1</sup>AB Sciex LP, 71 Four Valley Drive, Concord, Ontario L4K 4V8, Canada

<sup>2</sup>AB Sciex, 1201 Radio Road, Redwood City, CA 94065, USA

<sup>3</sup>Phenomenex Inc., 411 Madrid Ave., Torrance, CA 90501, USA

### Introduction

This technical note demonstrates the sensitive quantification of small molecule pharmaceutical compounds extracted from human plasma using minimal sample preparation and negative ion mode-based analysis. Lower limits of quantification (LLOQs) of 3.13 pg/mL and 25 pg/mL were achieved for Warfarin and Furosemide, respectively. Quantitative performance of the assay highlighted outstanding precision, accuracy, and linearity. Enhanced assay sensitivity was achieved through the application of a high-end triple quadrupole mass spectrometer to meet the demands of routine bioanalysis in complex matrices.

Demand for improved sensitivity in bioanalytical assays continues to increase as drug discovery and development programs focus on more efficacious, lower dosage compounds and as throughput demands drive the simplification of sample extraction and LC and MS methods. In many cases, the use of a more sensitive mass spectrometer is the easiest way to meet these needs. Having a system that offers technological improvements that deliver sensitivity gains across the mass range and in both polarities offers the bioanalytical scientist maximum flexibility to address the challenges outlined above. Here, Warfarin and Furosemide were extracted from human plasma and quantified using a negative ion mode approach on the SCIEX 7500 system (Figure 1). The improved front-end technology of the system enabled greater ion generation, capture and transmission to improve sensitivity for routine bioanalysis.

### Sample Preparation

Warfarin and Furosemide were spiked into 100 µL aliquots of human plasma at concentrations ranging from 3.13 to 100,000 pg/mL. Samples were extracted using protein precipitation with 300 µL Acetonitrile. The samples were then vortexed and centrifuged at 10,000 rpm for 25 minutes. The supernatant was collected for analysis.

### LC Conditions

**Column:** Kinetex™ 1.7 µm C18

**Dimensions:** 50 x 2.1 mm

**Part No.:** [00B-4475-AN](#)

**Mobile Phase:** A: 0.01 % Formic Acid in Water

B: 0.01 % Formic Acid in Acetonitrile

Gradient:	Time (min)	%B
	0	15
	0.25	15
	3	50
	3.1	95
	4	95
	4.1	15
	5	15

**Flow Rate:** 0.6 mL/min

**Injection Volume:** 1 µL

**Temperature:** Ambient

**LC System:** SCIEX® ExionLC™

**Detection:** MS/MS

**Detector:** SCIEX Triple Quad™ 7500

### MS/MS Conditions

**Polarity:** Negative

**Source Temperature:** 650 °C

**GS1:** 60 psi

**GS2:** 70 psi

**CUR:** 40 psi

**CAD:** 10

**IS:** 1700 V

### MS/MS Transitions and Parameters

Analyte	Q1 (m/z)	Q3 (m/z)	EP (V)	CE (V)	CXP (V)
Warfarin	307.1	250.1	-10	-31	-9
Furosemide	329	77.9	-10	-40	-9



Results and Discussion

Calibration curves were acquired across the concentration range of 3.13 to 100,000 pg/mL. Each concentration was analyzed in triplicate to assess method reproducibility. Strong linearity was observed across the concentration ranges analyzed, as demonstrated in **Figures 2 and 3** for Warfarin and Furosemide, respectively. **Table 1** summarizes the quantification results, including accuracy and precision. Excellent %CVs were achieved across all concentration levels with no interference in the blank human plasma samples for Warfarin and Furosemide. Using a generic sample preparation technique for the analysis of 100 µL of human plasma and a total run time of 5 minutes, the method provided LLOQs of 3.13 pg/mL and 25 pg/mL for Warfarin and Furosemide, respectively (**Figures 4 and 5**).

Figure 1. Extracted Ion Chromatograms (XICs) Representing the Matrix Blank and LLOQs of Warfarin and Furosemide Extracted from Human Plasma.

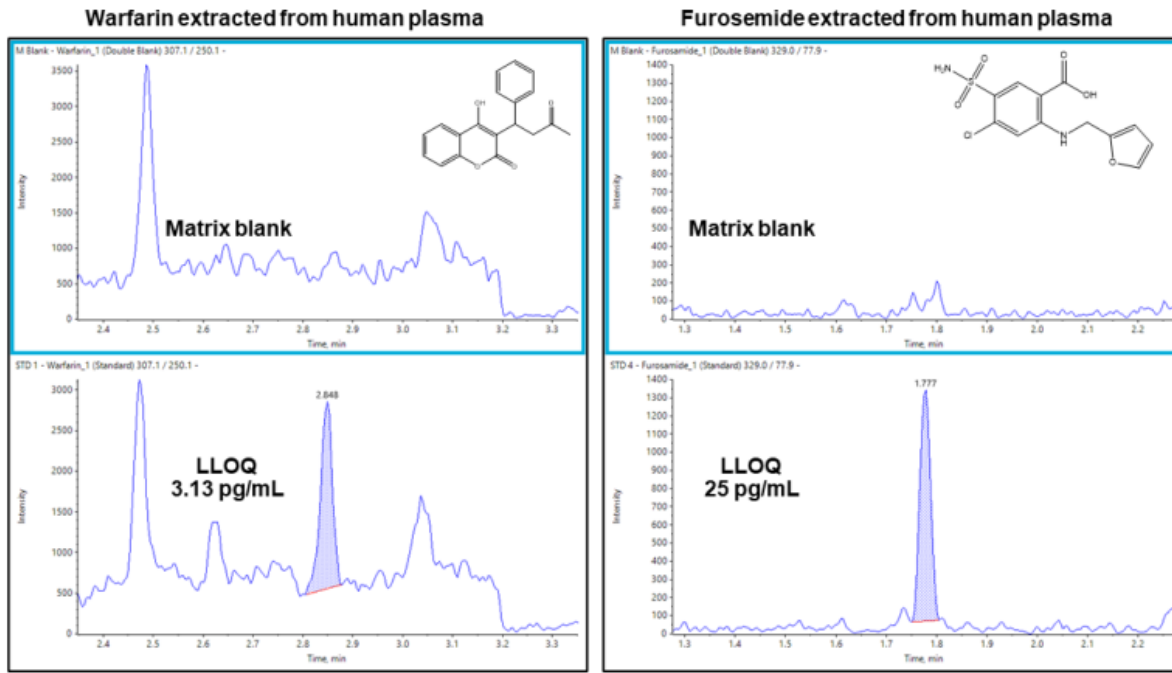


Figure 2. The Calibration Curve for Quantification of Warfarin in Human Plasma.


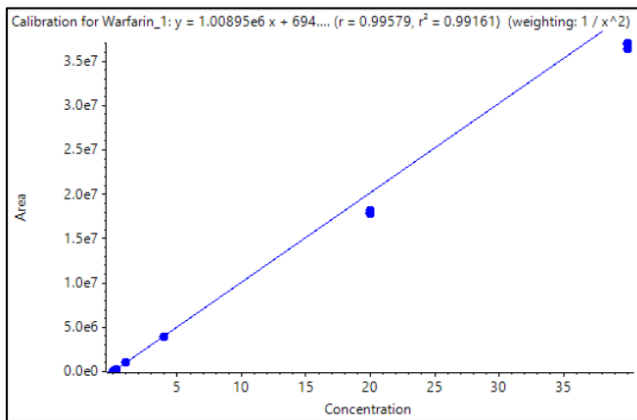



Figure 3. The Calibration Curve for the Quantification of Furosemide from Human Plasma.

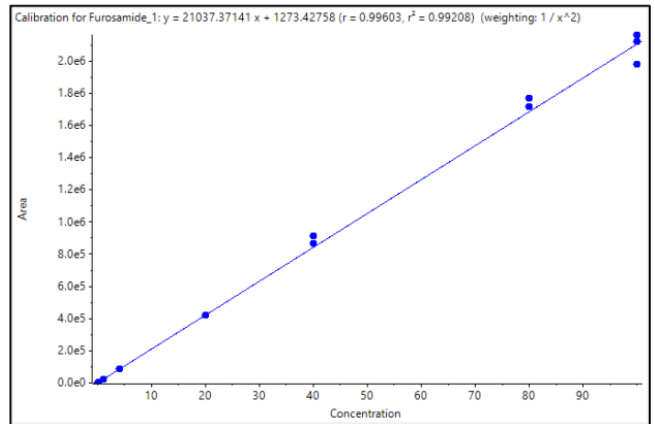


**Table 1.** Quantification Summary for Warfarin and Furosemide.

Concentration (fmol/ $\mu$ L)	Warfarin		Furosemide	
	Accuracy (%)	CV (%)	Accuracy (%)	CV (%)
3.13	100	9.2	N/A	N/A
6.25	94.5	4.1	N/A	N/A
12.5	103	1.4	N/A	N/A
25.0	106	2.6	104	15.3
62.5	115	1.7	91.6	10.2
250	103	0.8	88.5	1.6
1000	100	1.5	102	1.3
4000	97.1	1.4	104	1.2
20,000	89.1	1.3	99.4	0.1
40,000	91.2	0.9	107	3.1
80,000	N/A	N/A	104	1.8
100,000	N/A	N/A	99.2	4.5

**Figure 4.** XICs Demonstrating Warfarin Extraction from Human Plasma. a) Matrix Blank, and Warfarin present at b) 3.13 pg/mL, c) 6.25 pg/mL, d) 12.5 pg/mL, and e) 25 pg/mL.

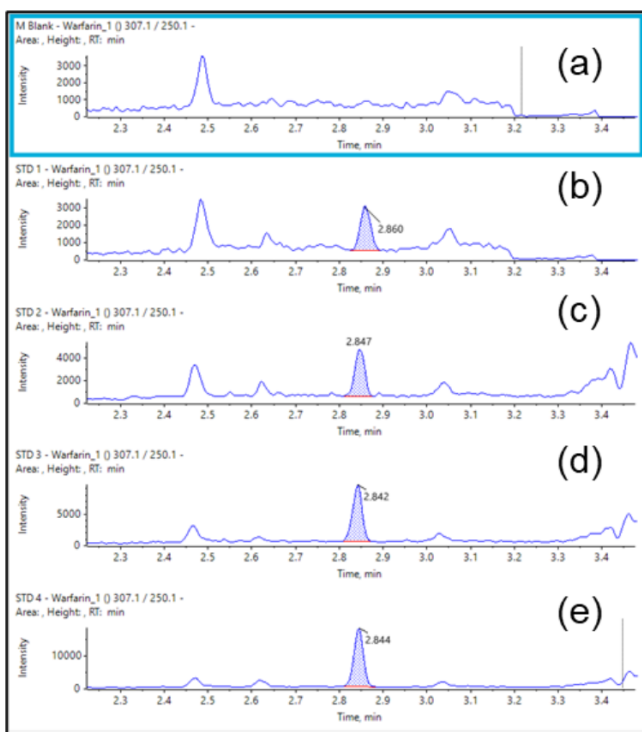

**Figure 5.** XICs Demonstrating Furosemide Extraction from Human Plasma. a) Matrix Blank, and Furosemide present at b) 25 pg/mL, c) 62.5 pg/mL, d) 250 pg/mL, and e) 1000 pg/mL.

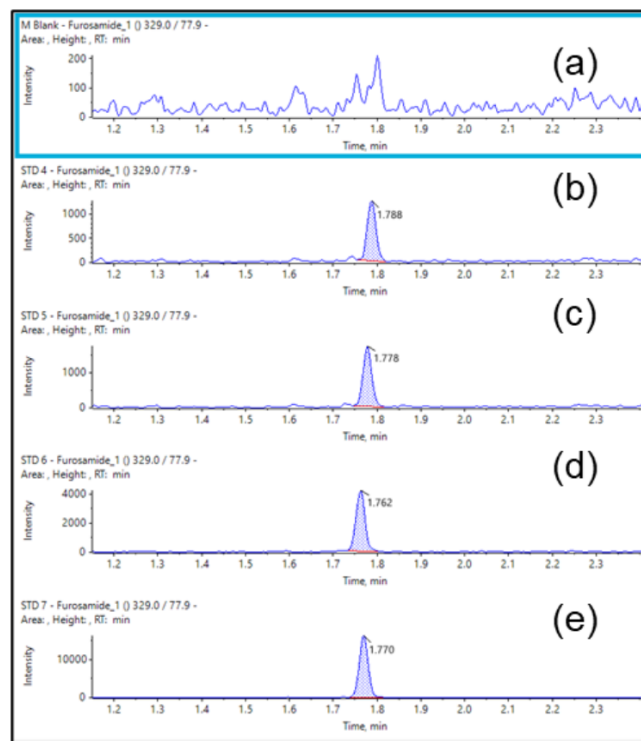

## Conclusions

Low-pg/mL level LLOQs were reached for Warfarin and Furosemide extracted from human plasma using minimal sample preparation. A highly sensitive assay for the quantification of small molecule pharmaceutical compounds requiring negative ion mode-based analysis was demonstrated on the SCIEX® 7500 system. The method demonstrated excellent accuracy, precision, and linearity at all concentration levels. A single platform for streamlined data acquisition, processing, and management with SCIEX OS software was presented. Overall, the SCIEX 7500 system enables pharmaceutical researchers the maximum flexibility to explore lower dosage, higher efficacy compounds, and improve the efficiency of bioanalysis.

## Kinetex™ Ordering Information

1.7 µm Minibore Columns (mm)	SecurityGuard™ ULTRA Cartridges†				
	30 x 2.1	50 x 2.1	100 x 2.1	150 x 2.1	3/pk
EVO C18	—	<a href="#">00B-4726-AN</a>	<a href="#">00D-4726-AN</a>	<a href="#">00F-4726-AN</a>	<a href="#">AJ0-9298</a>
Biphenyl	<a href="#">00A-4628-AN</a>	<a href="#">00B-4628-AN</a>	<a href="#">00D-4628-AN</a>	<a href="#">00F-4628-AN</a>	<a href="#">AJ0-9209</a>
XB-C18	<a href="#">00A-4498-AN</a>	<a href="#">00B-4498-AN</a>	<a href="#">00D-4498-AN</a>	<a href="#">00F-4498-AN</a>	<a href="#">AJ0-8782</a>
C18	<a href="#">00A-4475-AN</a>	<a href="#">00B-4475-AN</a>	<a href="#">00D-4475-AN</a>	<a href="#">00F-4475-AN</a>	<a href="#">AJ0-8782</a>
C8	<a href="#">00A-4499-AN</a>	<a href="#">00B-4499-AN</a>	<a href="#">00D-4499-AN</a>	<a href="#">00F-4499-AN</a>	<a href="#">AJ0-8784</a>
HILIC	<a href="#">00A-4474-AN</a>	<a href="#">00B-4474-AN</a>	<a href="#">00D-4474-AN</a>	—	<a href="#">AJ0-8786</a>
Phenyl-Hexyl	—	<a href="#">00B-4500-AN</a>	<a href="#">00D-4500-AN</a>	<a href="#">00F-4500-AN</a>	<a href="#">AJ0-8788</a>
F5	—	<a href="#">00B-4722-AN</a>	<a href="#">00D-4722-AN</a>	<a href="#">00F-4722-AN</a>	<a href="#">AJ0-9322</a>

for 2.1 mm ID

†SecurityGuard ULTRA Cartridges require holder, Part No.: [AJ0-9000](#)



**Need a different column size or sample preparation format?**

No problem! We have a majority of our available dimensions up on [www.phenomenex.com](http://www.phenomenex.com), but if you can't find what you need right away, our super helpful Technical Specialists can guide you to the solution via our online chat portal [www.phenomenex.com/Chat](http://www.phenomenex.com/Chat).

**Australia**

t: +61 (0)2-9428-6444  
auinfo@phenomenex.com

**Austria**

t: +43 (0)1-319-1301  
anfrage@phenomenex.com

**Belgium**

t: +32 (0)2 503 4015 (French)  
t: +32 (0)2 511 8666 (Dutch)  
beinfo@phenomenex.com

**Canada**

t: +1 (800) 543-3681  
info@phenomenex.com

**China**

t: +86 400-606-8099  
cninfo@phenomenex.com

**Czech Republic**

t: +420 272 017 077  
cz-info@phenomenex.com

**Denmark**

t: +45 4824 8048  
nordicinfo@phenomenex.com

**Finland**

t: +358 (0)9 4789 0063  
nordicinfo@phenomenex.com

**France**

t: +33 (0)1 30 09 21 10  
franceinfo@phenomenex.com

**Germany**

t: +49 (0)6021-58830-0  
anfrage@phenomenex.com

**Hong Kong**

t: +852 6012 8162  
hkinfo@phenomenex.com

**India**

t: +91 (0)40-3012 2400  
indiainfo@phenomenex.com

**Indonesia**

t: +62 21 5019 9707  
indoinfo@phenomenex.com

**Ireland**

t: +353 (0)1 247 5405  
eireinfo@phenomenex.com

**Italy**

t: +39 051 6327511  
italiainfo@phenomenex.com

**Japan**

t: +81 (0) 120-149-262  
jpinfo@phenomenex.com

**Luxembourg**

t: +31 (0)30-2418700  
nlinfo@phenomenex.com

**Mexico**

t: 01-800-844-5226  
tecnicomx@phenomenex.com

**The Netherlands**

t: +31 (0)30-2418700  
nlinfo@phenomenex.com

**New Zealand**

t: +64 (0)9-4780951  
nzinfo@phenomenex.com

**Norway**

t: +47 810 02 005  
nordicinfo@phenomenex.com

**Poland**

t: +48 22 104 21 72  
pl-info@phenomenex.com

**Portugal**

t: +351 221 450 488  
ptinfo@phenomenex.com

**Singapore**

t: +65 6559 4364  
sginfo@phenomenex.com

**Slovakia**

t: +420 272 017 077  
sk-info@phenomenex.com

**Spain**

t: +34 91-413-8613  
espinfo@phenomenex.com

**Sweden**

t: +46 (0)8 611 6950  
nordicinfo@phenomenex.com

**Switzerland**

t: +41 (0)61 692 20 20  
swissinfo@phenomenex.com

**Taiwan**

t: +886 (0) 0801-49-1246  
twinfo@phenomenex.com

**Thailand**

t: +66 (0) 2 566 0287  
thaiinfo@phenomenex.com

**United Kingdom**

t: +44 (0)1625-501367  
ukinfo@phenomenex.com

**USA**

t: +1 (310) 212-0555  
[www.phenomenex.com/chat](http://www.phenomenex.com/chat)

🌐 **All other countries/regions  
Corporate Office USA**

t: +1 (310) 212-0555  
[www.phenomenex.com/chat](http://www.phenomenex.com/chat)

**www.phenomenex.com**

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at [international@phenomenex.com](mailto:international@phenomenex.com)

**BE-HAPPY™  
GUARANTEE**

Your happiness is our mission. Take 45 days to try our products. If you are not happy, we'll make it right.

[www.phenomenex.com/behappy](http://www.phenomenex.com/behappy)

**Terms and Conditions**

Subject to Phenomenex Standard Terms and Conditions, which may be viewed at [www.phenomenex.com/phx-terms-and-conditions-of-sale](http://www.phenomenex.com/phx-terms-and-conditions-of-sale).

**Trademarks**

Kinetex, SecurityGuard, and BE-HAPPY are trademarks of Phenomenex. SCIEIX is a registered trademark and Triple Quad and ExionLC are trademarks of AB SCIEIX Pte. Ltd.

**Disclaimer**

Comparative separations may not be representative of all applications.

SecurityGuard is patented by Phenomenex. U.S. Patent No. 6,162,362

**CAUTION: this patent only applies to the analytical-sized guard cartridge holder, and does not apply to SemiPrep, PREP, or ULTRA holders, or to any cartridges.**

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures.

© 2022 Phenomenex, Inc. All rights reserved.

