

## Separation of Miconazole Nitrate and its Organic Impurities per USP Monograph using Kinetex™ 2.6 µm Phenyl-Hexyl Column

Lauren Nakasone, Bryan Tackett, PhD, and Zeshan Aqeel  
Phenomenex Inc., 411 Madrid Ave., Torrance, CA 90501, USA

### Overview

Miconazole Nitrate belongs to the imidazole group of antifungal agents. The development of a quick and efficient analysis of Miconazole Nitrate and its related organic impurities is of interest for drug manufacturers. In this application note, we report the separation of Miconazole Nitrate and its related organic impurities using the Kinetex 2.6 µm Phenyl-Hexyl column specified by the USP monograph for Miconazole Nitrate.

System suitability per USP Monograph for the Miconazole Nitrate Assay is a resolution no less than (NLT) 1.5 between Miconazole Nitrate and Miconazole Related Compound F, a symmetry factor no more than (NMT) 2.0 for Miconazole Nitrate, and a percent relative standard deviation (%RSD) of NMT 0.73 % for Miconazole Nitrate. All system suitability requirements for Miconazole Nitrate Assay were met using the Kinetex Phenyl-Hexyl column (Figure 1 and 2).

System suitability per USP Monograph for the Miconazole Nitrate Organic Impurities is a resolution NLT 1.5 between Miconazole Related Compound C and Miconazole Related Compound I, a resolution NLT 1.5 between Miconazole Related Compound I and Econazole Nitrate, a resolution NLT 1.5 between Miconazole Related Compound F and Miconazole Nitrate, as well as a %RSD NMT 3.0 % for Miconazole Nitrate. The separation of Miconazole Nitrate and its organic impurities was achieved with all requirements for System Suitability for Organic Impurities met using the Kinetex Phenyl-Hexyl column (Figure 3).

All solutions were prepared as indicated in the USP Monograph for Miconazole Nitrate. USP Miconazole Nitrate RS (Catalog No. 1443500), USP Econazole Nitrate RS (Catalog No. 1231808), USP Miconazole Related Compound C RS (Catalog No. 1443431), USP Miconazole Related Compound F RS (Catalog No. 1443464), and USP Miconazole Related Compound I RS (Catalog No. 1443497) were purchased from USP.

### LC-UV Conditions

**Column:** Kinetex 2.6 µm Phenyl-Hexyl

**Dimension:** 100 x 4.6 mm

**Part No.:** [00D-4495-E0](#)

**Mobile Phase:** A: Methanol/Water/1 M Triethylammonium Acetate (30:70:1, v/v/v)

B: Acetonitrile/Methanol/1 M Triethylammonium Acetate (25:75:0.2, v/v/v)

Gradient:	Time (min)	%B
	0	30
	5	30
	10	56
	27	56
	30	75
	35	75
	36	30
	40	30

**Flow Rate:** 0.8 mL/min

**Injection Volume:** 10 µL

**Temperature:** 40 °C

**Detector:** UV @ 215


**System:** Waters® ACQUITY® I-Class UPLC

**Table 1.** Preparation of Solutions

Solution	Composition
Diluent – Assay	Methanol/Water (70:30, v/v)
Standard Solution – Assay	0.1 mg/mL USP Miconazole Nitrate RS in Diluent
System Suitability Solution – Assay	0.1 mg/mL of USP Miconazole Nitrate RS and 6 µg/mL of USP Miconazole Related Compound F RS in Diluent
Standard Solution – Organic Impurities	1.2 µg/mL each of USP Miconazole Nitrate RS, USP Econazole Nitrate RS, USP Miconazole Related Compound C RS, USP Miconazole Related Compound F RS, and USP Miconazole Related Compound I RS in Diluent




**Figure 1. Standard Solution – Assay**

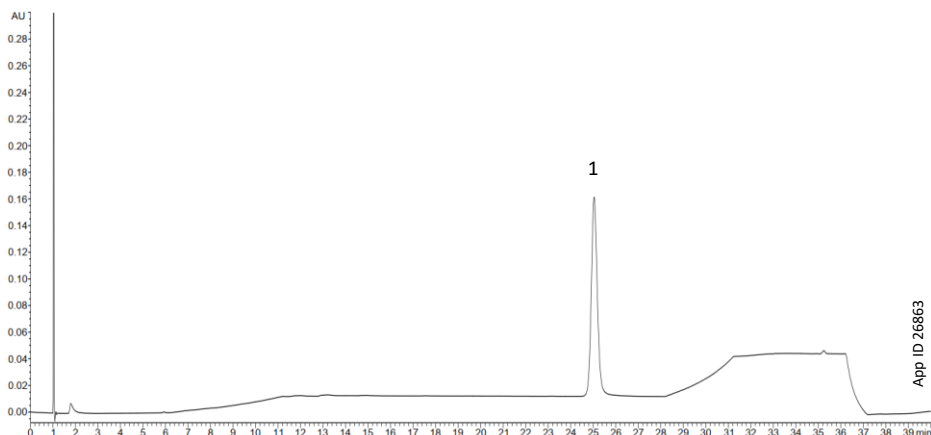

Peak No.	Analyte	Symmetry Factor	Area %RSD
1	Miconazole Nitrate	1.11	0.421
N=6 Injections			

**Figure 2. System Suitability Solution – Assay**

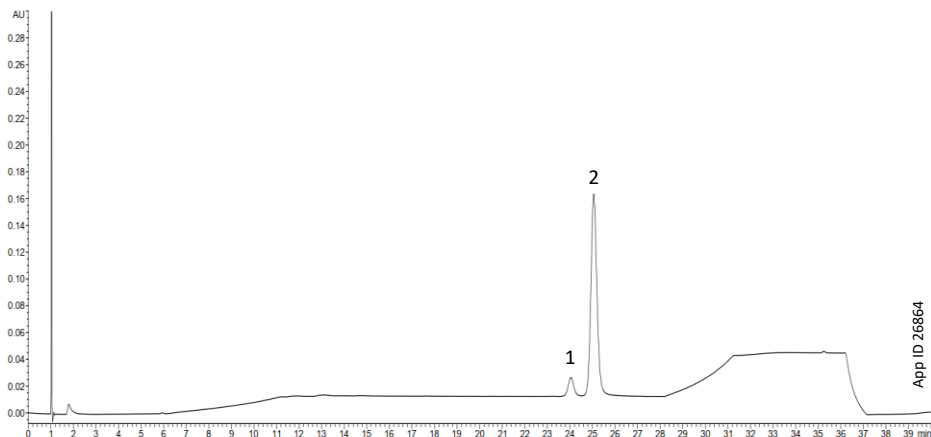

Peak No.	Analyte	Resolution
1	Miconazole Related Compound F	2.1
2	Miconazole Nitrate	
N=6 Injections		



**Figure 3.** Standard Solution – Organic Impurities



Peak No.	Analyte	Resolution	Area %RSD
1	Miconazole Related Compound C	-	-
2	Miconazole Related Compound I	2.417	-
3	Econazole Nitrate	-	-
4	Miconazole Related Compound F	-	-
5	Miconazole Nitrate	2.2	2.592

N=6 Injections



**Need a different column size or sample preparation format?**

No problem! We have a majority of our available dimensions up on [www.phenomenex.com](http://www.phenomenex.com), but if you can't find what you need right away, our super helpful Technical Specialists can guide you to the solution via our online chat portal [www.phenomenex.com/LiveChat](http://www.phenomenex.com/LiveChat).

**Australia**

t: +61 (0)2-9428-6444  
auiinfo@phenomenex.com

**Austria**

t: +43 (0)1-319-1301  
anfrage@phenomenex.com

**Belgium**

t: +32 (0)2 503 4015 (French)  
t: +32 (0)2 511 8666 (Dutch)  
beinfo@phenomenex.com

**Canada**

t: +1 (800) 543-3681  
info@phenomenex.com

**China**

t: +86 400-606-8099  
cninfo@phenomenex.com

**Czech Republic**

t: +420 272 017 077  
cz-info@phenomenex.com

**Denmark**

t: +45 4824 8048  
nordicinfo@phenomenex.com

**Finland**

t: +358 (0)9 4789 0063  
nordicinfo@phenomenex.com

**France**

t: +33 (0)1 30 09 21 10  
franceinfo@phenomenex.com

**Germany**

t: +49 (0)6021-58830-0  
anfrage@phenomenex.com

**Hong Kong**

t: +852 6012 8162  
hkinfo@phenomenex.com

**India**

t: +91 (0)40-3012 2400  
indiainfo@phenomenex.com

**Indonesia**

t: +62 21 5010 9707  
indoinfo@phenomenex.com

**Ireland**

t: +353 (0)1 247 5405  
eireinfo@phenomenex.com

**Italy**

t: +39 051 6327511  
italiainfo@phenomenex.com

**Japan**

t: +81 (0) 120-149-262  
jpinfo@phenomenex.com

**Luxembourg**

t: +31 (0)30-2418700  
nlinfo@phenomenex.com

**Mexico**

t: 01-800-844-5226  
tecnicomx@phenomenex.com

**The Netherlands**

t: +31 (0)30-2418700  
nlinfo@phenomenex.com

**New Zealand**

t: +64 (0)9-4780951  
nzinfo@phenomenex.com

**Norway**

t: +47 810 02 005  
nordicinfo@phenomenex.com

**Poland**

t: +48 22 104 21 72  
pl-info@phenomenex.com

**Portugal**

t: +351 221 450 488  
ptinfo@phenomenex.com

**Singapore**

t: +65 800-852-3944  
sginfo@phenomenex.com

**Slovakia**

t: +420 272 017 077  
sk-info@phenomenex.com

**Spain**

t: +34 91-413-8613  
espinfo@phenomenex.com

**Sweden**

t: +46 (0)8 611 6950  
nordicinfo@phenomenex.com

**Switzerland**

t: +41 (0)61 692 20 20  
swissinfo@phenomenex.com

**Taiwan**

t: +886 (0) 0801-49-1246  
twinfo@phenomenex.com

**Thailand**

t: +66 (0) 2 566 0287  
thaiinfo@phenomenex.com

**United Kingdom**

t: +44 (0)1625-501367  
ukinfo@phenomenex.com

**USA**

t: +1 (310) 212-0555  
info@phenomenex.com

🌐 **All other countries/regions  
Corporate Office USA**

t: +1 (310) 212-0555  
info@phenomenex.com

**www.phenomenex.com**

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at [international@phenomenex.com](mailto:international@phenomenex.com)

**BE-HAPPY™  
GUARANTEE**

Your happiness is our mission. Take 45 days to try our products. If you are not happy, we'll make it right.

[www.phenomenex.com/behappy](http://www.phenomenex.com/behappy)

**Terms and Conditions**

Subject to Phenomenex Standard Terms and Conditions, which may be viewed at [www.phenomenex.com/TermsAndConditions](http://www.phenomenex.com/TermsAndConditions).

**Trademarks**

Kinetex and BE-HAPPY are trademarks of Phenomenex. Waters and ACQUITY are registered trademarks of Waters Technologies Corporation.

**Disclaimer**

Phenomenex is in no way affiliated with Waters Technologies Corporation.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures.

© 2022 Phenomenex, Inc. All rights reserved.

