

Intact Identification of Adeno-Associated Virus Capsid Proteins

Lorne Nelson, Mastroeh Chamanian, PhD, Roxana Eggleston-Rangel, and Bryan Tackett, PhD
Phenomenex, Inc., 411 Madrid Ave., Torrance, CA 90501 USA

Overview

The use of Adeno-Associated Viruses (AAV) as an effective vector for the delivery of ssDNA in gene therapy continues to gain momentum. A chief attribute to this success is the broad tissue tropism enabled by a range of serotypes. To date, approximately 13 different serotypes have been identified with a range of shared amino acid sequence identity. Structurally, the viral capsid is a sixty-protein icosahedron comprised of three capsid, or viral proteins (VP), non-covalently assembled in a ratio of 1:1:10 (VP1:VP2:VP3).

These three proteins share a high degree of sequence homology, as VP2 and VP3 represent iterative, N-terminal truncations of the VP1 protein. The resulting approximate masses are VP1 – 81 kDa, VP2 – 66 kDa, and VP3 – 60 kDa. Given the inter-serotype and structural protein sequence similarities, protein identification, relative protein ratio, and post-translational modification assessments are critical. However, as the complete capsid can range in mass from 3.67 MDa, for an empty capsid, to 4.75 MDa for a capsid containing its payload (up to 4.7kb ssDNA), capsid denaturation and dissociation are required prior to interrogation.

In this application note, we profile denatured and dissociated capsid proteins from AAV5 and AAV9, via analysis with LC-MS using a Biozen™ Intact XB-C8 reversed phase column. It is notable that, although no covalent attachments are present within the capsid structure, denaturation without reduction can lead to a covalent complex of VP1 and VP2 in AAV9.

Sample Preparation

Step	Description
Sample:	AAV with stock concentration of 2E13 vg/mL (approx. 0.122 µg/µL protein)
Dilution:	5X dilution with 1X Dulbecco's Phosphate Buffered Saline
Denaturation:	10 % v/v 100 mM Acetic or Formic Acid (as indicated)
Reduction:	500 mM Tris (2-carboxyethyl) Phosphine for a 30 mM final concentration (as indicated)
Incubation:	22 °C for 15 min
Centrifugation:	9000 rpm for 5 min
Transfer:	Transfer to HPLC sample vial

LC Conditions

Column:	Biozen 3.6 µm Intact XB-C8	
Dimension:	150 x 2.1 mm	
Part No.:	<u>00F-4766-AN</u>	
Mobile Phase:	A: 0.1 % Trifluoroacetic Acid in Water B: 0.1 % Trifluoroacetic Acid in Acetonitrile	
Gradient:	Time (min)	%B
	0	10
	3	10
	3.1	35
	7.1	35
	21.94	65
	22.04	90
	25.04	90
	25.14	10
	29.14	10
Flow Rate:	0.3 mL/min	
Injection Volume:	5 µL	
Temperature:	60 °C	
Instrument:	Vanquish™ UHPLC	
Detection:	MS	
Detector:	Q Exactive™ Orbitrap™ Plus	

MS Conditions

Scan Type:	Full MS
Resolution:	70,000
AGC Target:	3e6 ms
Maximum IT:	100 ms
Scan Range:	500 to 3500 m/z



Figure 1. Total Ion Chromatogram of AAV5 Capsid Proteins Denatured with Acetic Acid, at a Concentration of 3.2E12 vg/mL (Approx. 98 ng Protein).

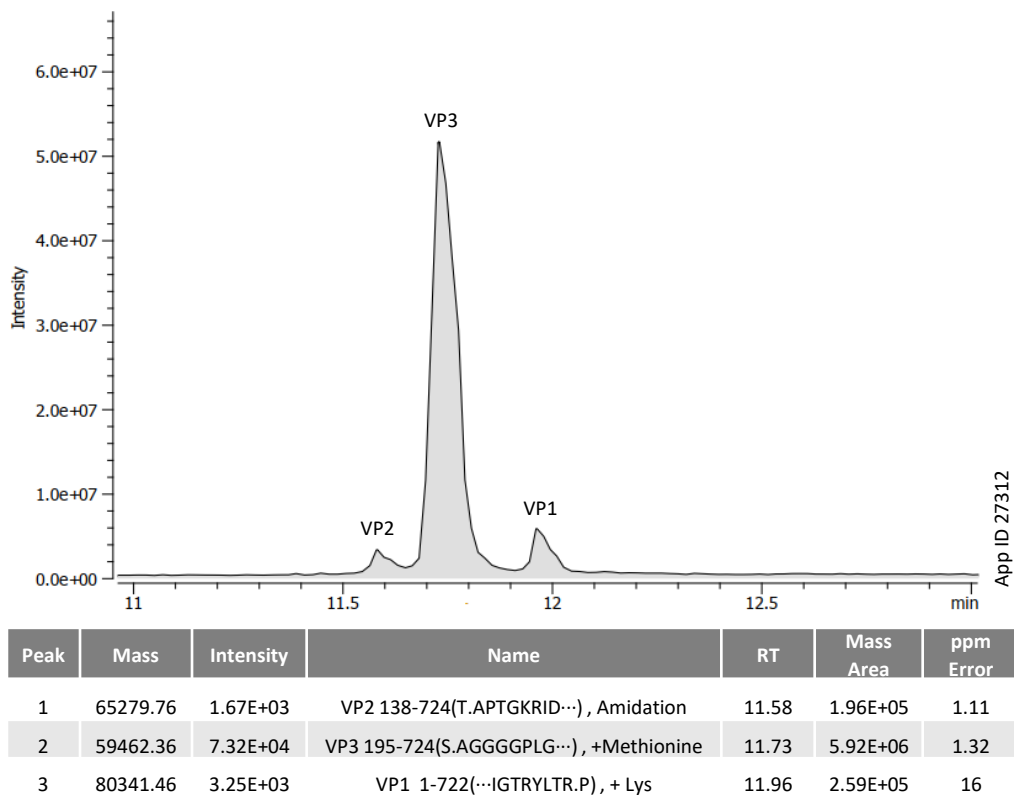


Figure 2. Total Ion Chromatogram of AAV9 Capsid Proteins Denatured with Acetic Acid and Reduced with Tris (2-carboxyethyl) Phosphine, at a Concentration of 3.2E12 vg/mL (Approx. 92 ng Protein).


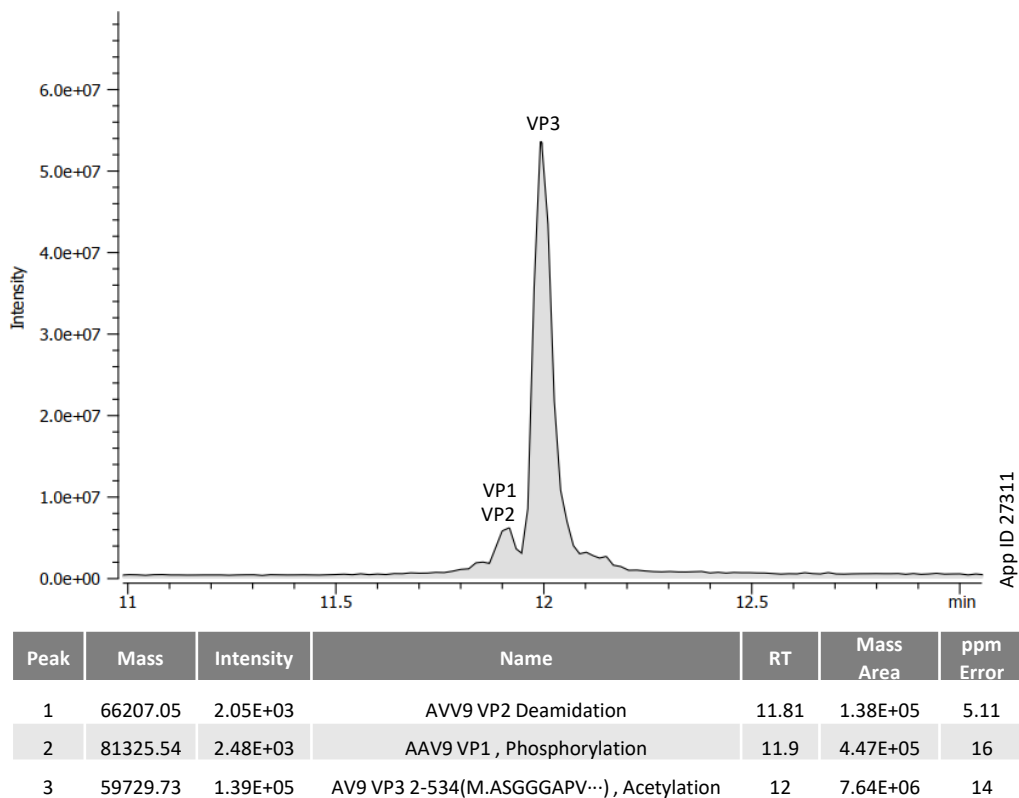


Figure 3. Total Ion Chromatogram of AAV9 Capsid Proteins Denatured with Formic Acid and Reduced with Tris (2-carboxyethyl) Phosphine, at a Concentration of 3.2E12 vg/mL (Approx. 92 ng Protein).



Peak	Mass	Intensity	Name	RT	Mass Area	ppm Error
1	66207.05	9.58E+03	AV9 VP2 , Deamidation	11.84	5.23E+05	4.99
2	81279.45	8.27E+03	AAV9 VP1 1-733(...IGTRYLTR.N) , +2Met	11.92	1.19E+06	4.43
3	59685.96	5.43E+04	AAV9 VP3 203-735(M.ASGGGAPV...), Amidation	11.99	1.10E+07	6.54

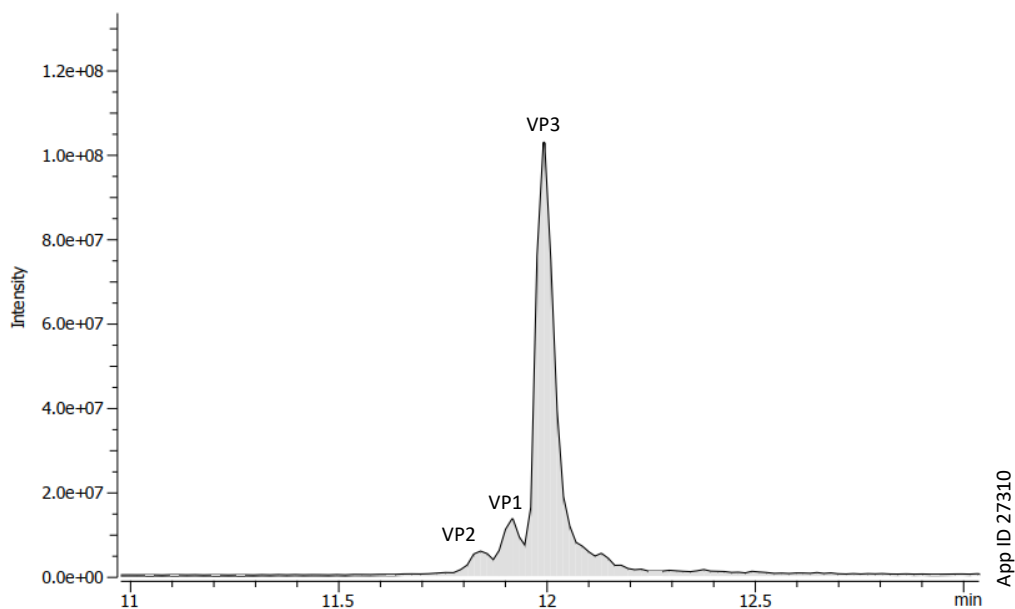

Need a different column size or sample preparation format?

No problem! We have a majority of our available dimensions up on www.phenomenex.com, but if you can't find what you need right away, our super helpful Technical Specialists can guide you to the solution via our online chat portal www.phenomenex.com/Chat.

Australia

t: +61 (0)2-9428-6444
auinfo@phenomenex.com

Austria

t: +43 (0)1-319-1301
anfrage@phenomenex.com

Belgium

t: +32 (0)2 503 4015 (French)
t: +32 (0)2 511 8666 (Dutch)
beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681
info@phenomenex.com

China

t: +86 400-606-8099
cninfo@phenomenex.com

Czech Republic

t: +420 272 017 077
cz-info@phenomenex.com

Denmark

t: +45 4824 8048
nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063
nordicinfo@phenomenex.com

France

t: +33 (0)1 30 09 21 10
franceinfo@phenomenex.com

Germany

t: +49 (0)6021-58830-0
anfrage@phenomenex.com

Hong Kong

t: +852 6012 8162
hkinfo@phenomenex.com

India

t: +91 (0)40-3012 2400
indiainfo@phenomenex.com

Indonesia

t: +62 21 5019 9707
indoinfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405
eireinfo@phenomenex.com

Italy

t: +39 051 6327511
italiainfo@phenomenex.com

Japan

t: +81 (0) 120-149-262
jpinfo@phenomenex.com

Luxembourg

t: +31 (0)30-2418700
nlinfo@phenomenex.com

Mexico

t: 01-800-844-5226
tecnicomx@phenomenex.com

The Netherlands

t: +31 (0)30-2418700
nlinfo@phenomenex.com

New Zealand

t: +64 (0)9-4780951
nzinfo@phenomenex.com

Norway

t: +47 810 02 005
nordicinfo@phenomenex.com

Poland

t: +48 22 104 21 72
pl-info@phenomenex.com

Portugal

t: +351 221 450 488
ptinfo@phenomenex.com

Singapore

t: +65 6559 4364
sginfo@phenomenex.com

Slovakia

t: +420 272 017 077
sk-info@phenomenex.com

Spain

t: +34 91-413-8613
espinfo@phenomenex.com

Sweden

t: +46 (0)8 611 6950
nordicinfo@phenomenex.com

Switzerland

t: +41 (0)61 692 20 20
swissinfo@phenomenex.com

Taiwan

t: +886 (0) 0801-49-1246
twinfo@phenomenex.com

Thailand

t: +66 (0) 2 566 0287
thaiinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367
ukinfo@phenomenex.com

USA

t: +1 (310) 212-0555
www.phenomenex.com/chat

**🌐 All other countries/regions
Corporate Office USA**

t: +1 (310) 212-0555
www.phenomenex.com/chat

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at international@phenomenex.com

BE-HAPPY™
GUARANTEE

Your happiness is our mission. Take 45 days to try our products. If you are not happy, we'll make it right.

www.phenomenex.com/behappy

Terms and Conditions

Subject to Phenomenex Standard Terms and Conditions, which may be viewed at www.phenomenex.com/phx-terms-and-conditions-of-sale.

Trademarks

Biozen and BE-HAPPY are trademarks of Phenomenex. Vanquish, Q Exactive, and Orbitrap are trademarks of Thermo Fisher Scientific, Inc.

Disclaimer

Comparative separations may not be representative of all applications. Phenomenex is in no way affiliated with Thermo Fisher Scientific, Inc.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures.

© 2023 Phenomenex, Inc. All rights reserved.

