



TN-1310

Micro-flow Peptide Separations

Roxana Eggleston-Rangel, Namrata Saxena, Jason Anspach, PhD, and Bryan Tackett, PhD
Phenomenex, Inc., 411 Madrid Ave., Torrance, CA 90501 USA

Introduction

Proteomics is the study of all proteins in an organism (proteome) or part of an organism. To study the proteome, Liquid Chromatography (LC) coupled to mass spectrometry (MS) has become the gold standard. Here, the role of LC is to provide the necessary resolution among peptides within a sample in order to be quantified or identified by MS.

An organism's proteome can be very complex. For example, a human cell contains about 10,000 protein producing coding genes, which in turn result in thousands of proteins with a myriad of different chemistries depending on their amino acid sequences and Post Translational Modifications (PTMs). In bottom-up proteomics, where an enzyme is used to digest the proteome prior to LC-MS analysis, understanding the chemistry of the sample being analyzed can be very beneficial. In this way, one can better choose the proper LC column phase chemistry to obtain the best resolution that provides the highest MS peptide identification and/or quantitation.

In this technical note, a mixture of 20 synthetic peptides with a wide mass and retention time range coverage (Table 1) was used to screen four micro-flow phase chemistries and gain insights into their peptide selectivity. A polar column that gives increased polar peptide retention, a biphenyl phase that provides aromatic selectivity, an F5 that gives aromatic plus electrostatic retention, and an XB-C18 with isobutyl side chains to improve peak shapes were used and behaved as expected.

Table 1. SCIEX PepCalMix Peptide Amino Acid Sequences Used and their Respective Monitored M+H Species.

Peptide Sequence	Analyte - m/z
AETSELHTSLK	408.5501+++
GAYVEVTAK	473.2602++
IGNEQGVSR	485.2530++
LDSTSIPVAK	519.7997++
AGLIVAEGVTK	533.3233++
LVGTPAEER	491.2656++
LGLDFDSFR	540.2734++
GFTAYYIPR	549.2863++
SGGLLWQLVR	569.8340++
AVGANPEQLTR	583.3136++
SAEGLDASASLR	593.8005++
VFTPLEVDVAK	613.3495++
VGNEIQYVALR	636.3527++
YIELAPGVDNSK	657.3450++
DGTFAVDGPVIAK	677.8583++
YDSINNTEVSGIR	739.3615++
SPYVITGPGVVEYK	758.9105++
ALENDIGVPSDATVK	768.9034++
AVFYAPQIPLYANK	589.3183+++
TVESLFPPEEAETPGSAVR	643.6540+++

LC Conditions

- Column:** Kinetex™ 2.6 µm Biphenyl
Luna™ Omega 3 µm Polar C18
Kinetex 2.6 µm XB-C18
Kinetex 2.6 µm F5
- Dimension:** 150 x 0.3 mm
- Part No.:** [00F-4622-AC](#) (Biphenyl)
[00F-4760-AC](#) (Polar C18)
[00F-4496-AC](#) (XB-C18)
[00F-4723-AC](#) (F5)
- Mobile Phase:** A: 0.1 % Formic Acid in Water
B: 0.1 % Formic Acid in Acetonitrile
- Gradient:**

Time (min)	%B
0	2
60	40
60.1	80
65	80
65.1	2
70	2
- Flow Rate:** 10 µL/min
- Injection:** 1 µL
- Temperature:** 40 °C
- Detector:** Q Exactive™ Plus Orbitrap™
- System:** NanoLC™ 425 (SCIEX®)
- Detection:** MS

MS Conditions

- Scan Type:** Full MS SIM
- Polarity:** Positive
- Resolution:** 70,000
- Scan Range:** 400 to 1650 m/z
- AGC Target:** 3e6
- Maximum IT:** 100 ms



Results and Discussion

A 20 isotopically labeled peptide mix covering a wide range of retention times and masses was used to screen 4 different micro column chemistry phases (Table 1). First, a Kinetex™ 2.6 µm XB-C18 stationary phase, that employs core-shell technology to reduce peptide diffusion and silanol groups present on the stationary phase, gave narrower peaks and higher peak capacities than the rest of the screened columns (Figure 1 and Table 2).

However, the first eluting peptide (IGNEQGVSR), which contains acidic groups like Glutamic Acid, was retained longer on a column with more polar character (Figure 2). In the case of peptides with aromatic character, the Kinetex Biphenyl column provided better resolution than the rest of the screened selectivities (Figures 3-6).

For example, when comparing AGLIVAEGVTK and YIELAPGVDNSK (Figures 3 and 4), YIELAPGVDNSK (with its aromatic Tyrosine) was fully resolved using a Kinetex Biphenyl column as opposed to the Kinetex XB-C18 or other selectivities. The same can be said of SGGLLWQLVR and AVYFYAPQIPLYANK (Figures 5 and 6). AVYFYAPQIPLYANK contains two Tyrosine and one Phenylalanine, making it more aromatic than the SGGLLWQLVR peptide. The ability to resolve aromatic rich peptides may prove useful in proteomics applications that deal with W,Y, F, and H rich proteins. Examples include breast cancer biomarkers, DNA/RNA-protein binding, carbohydrate-protein binding, and cation-n protein interactions.

Amino acid composition can vary widely and choosing a column can be more complicated than shown in the above examples. Therefore, testing more than one phase can be beneficial. For instance, in the case of the separation of GFTAYYIPR and DGTFAVDGPGVIAK (Figures 7 and 8), the Kinetex Biphenyl phase did not provide the best resolving power. The Kinetex F5 column provided the resolution needed to separate the more polar, charged, and aromatic peptide DGTFAVDGPGVIAK from GFTAYYIPR.

Figure 1. Extracted Ion Chromatogram (XIC) of PepCalMix Peptides on a Kinetex 2.6 µm XB-C18 column.

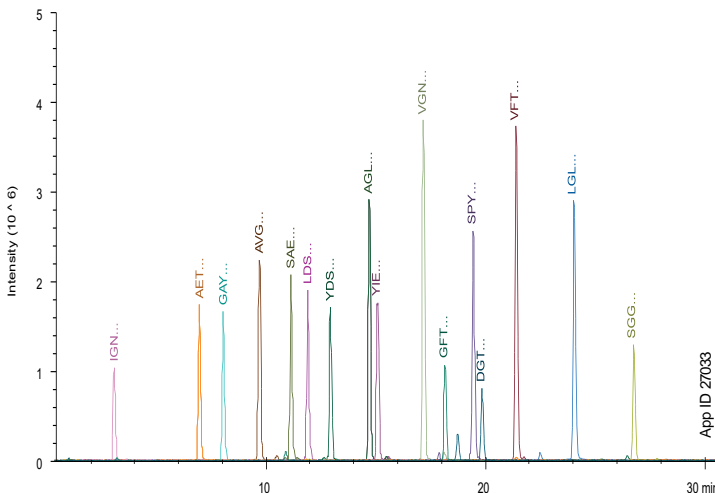


Table 2. Average Peak Widths and Peak Capacities for the 3 Kinetex Phases Used.

Column	Avg Peak Width	Peak Capacity
Kinetex 2.6 µm XB-C18	0.35	72
Kinetex 2.6 µm Biphenyl	0.42	68
Kinetex 2.6 µm F5	0.45	59

Figure 2. Retention Time Column Comparison Showing Retention Time Differences of Peptides when Using a Kinetex F5, a Kinetex Biphenyl, a Kinetex XB-C18 or a Luna™ Omega Polar Column.

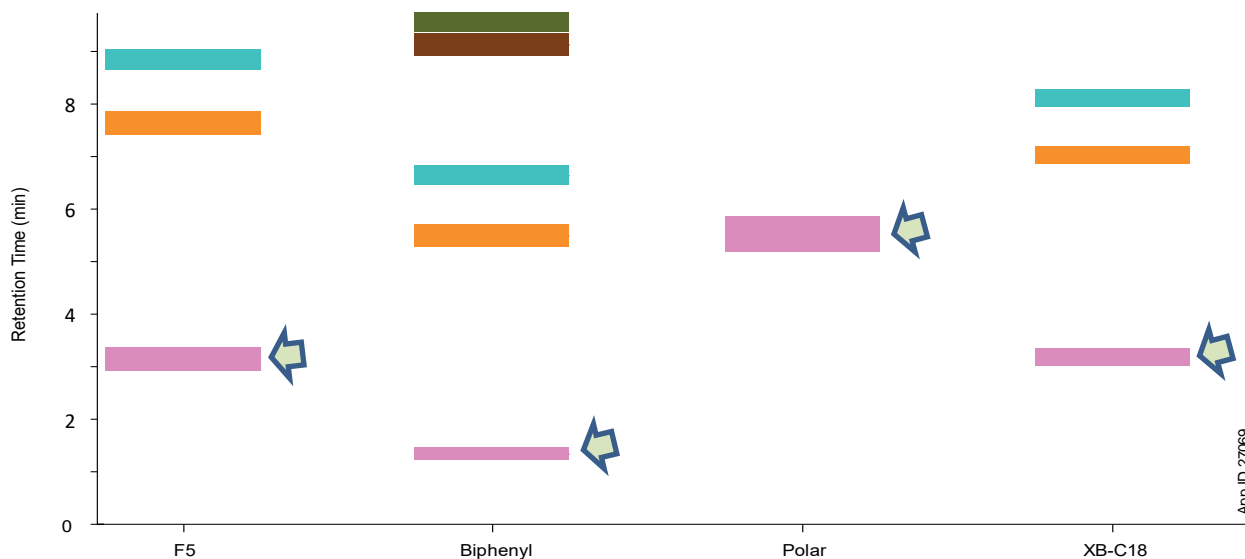


Figure 3. XIC of Peptides 1)AGLIVAEGVTK and 2)YIELAPGVDNSK on Kinetex™ 2.6 µm XB-C18 Column.

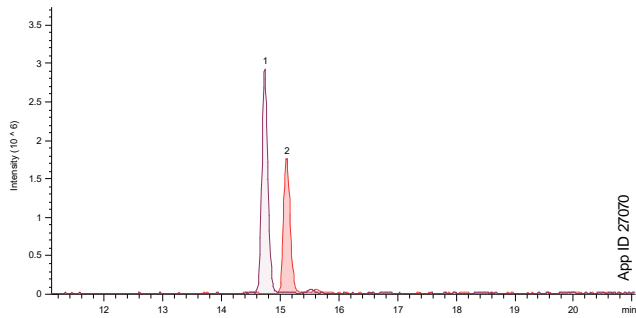


Figure 4. XIC of Peptides 1)AGLIVAEGVTK and 2)YIELAPGVDNSK on a Kinetex 2.6 µm Biphenyl Column.

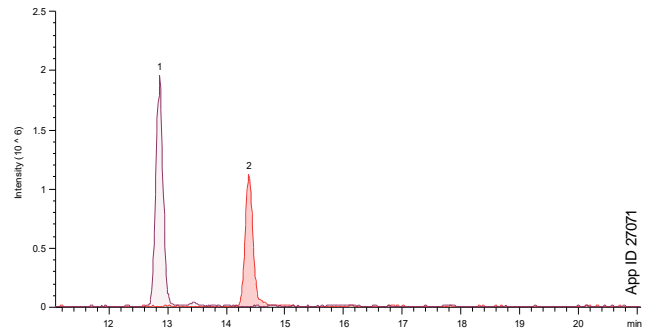


Figure 5. XIC of Peptides 1)SGLLWQLVR and 2)AVYFYAPQIPLYANK on Kinetex 2.6 µm XB-C18 Column.

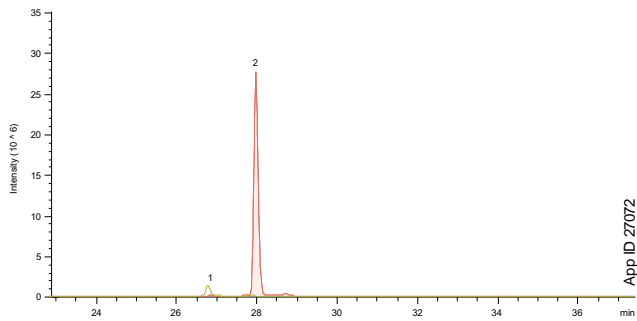


Figure 6. XIC of Peptides 1)SGLLWQLVR and 2)AVYFYAPQIPLYANK on Kinetex 2.6 µm Biphenyl Column.

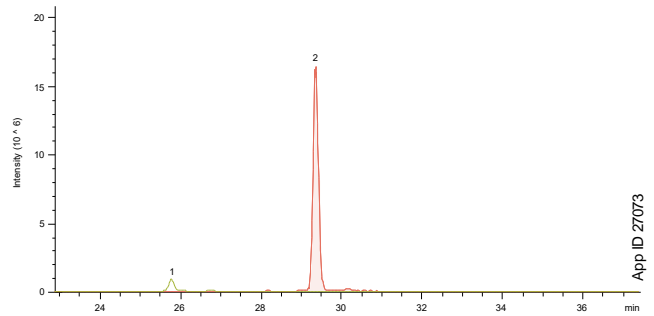


Figure 7. XIC of Peptides 1)GFTAYIPR and 2)DGTFAVDGPGVIAK on Kinetex 2.6 µm Biphenyl Column.


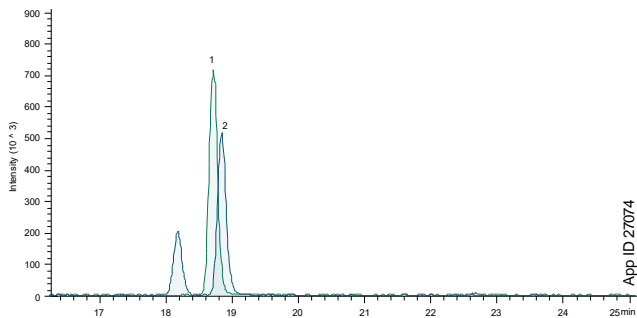


Figure 8. XIC of Peptides 1)DGTFAVDGPGVIAK and 2)GFTAYIPR on Kinetex 2.6 µm F5 Column.



Conclusions

Proteomics samples are composed of thousands of proteins, and in the case of bottom-up proteomics, thousands of peptides. These peptides are made of many different amino acid sequences in addition to PTMs that give those peptides unique chemistries. In order to get the most out of an LC-MS workflow, the column phase chemistry should be carefully explored. As shown here, when dealing with early eluters, a polar column may be the best choice. In the case of aromatic peptides, a Biphenyl or an F5 column could provide the optimal resolution. When in doubt, an XB-C18 column can be a good place to start to obtain optimal peak shape and determine specific needs depending on the sample chemistry.

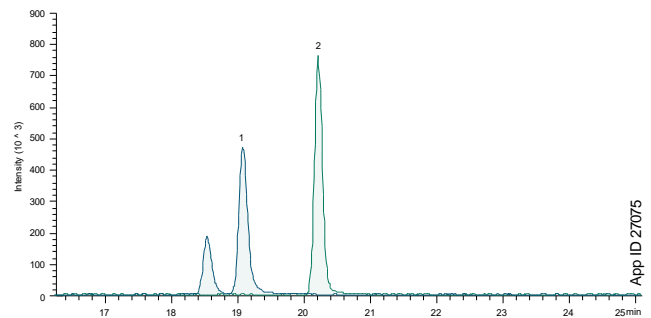

Kinetex™ Micro LC Columns Ordering Information

2.6 µm Micro LC Columns (mm)						
Part	Phases	30 x 0.3	50 x 0.3	100 x 0.3	150 x 0.3	50 x 0.5
XB-C18	00A-4496-AC	00B-4496-AC	00D-4496-AC	00F-4496-AC	00B-4496-AF	00F-4496-AF
Biphenyl	—	00B-4622-AC	—	00F-4622-AC	00B-4622-AF	—
C18	00A-4462-AC	00B-4462-AC	—	00F-4462-AC	00B-4462-AF	—
EVO C18	—	00B-4725-AC	—	00F-4725-AC	00B-4725-AF	—
F5	—	00B-4723-AC	00D-4723-AC	00F-4723-AC	00B-4723-AF	—

Luna™ Omega Micro LC Columns Ordering Information

3 µm Micro LC Columns (mm)							Trap Column
Phases	50 x 0.30	100 x 0.30	150 x 0.30	50 x 0.50	100 x 0.50	150 x 0.50	20 x 0.3
Polar C18	00B-4760-AC	00D-4760-AC	00F-4760-AC	00B-4760-AF	00D-4760-AF	00F-4760-AF	—
PS C18	00B-4758-AC	00D-4758-AC	00F-4758-AC	00B-4758-AF	00D-4758-AF	00F-4758-AF	05M-4758-AC



Need a different column size or sample preparation format?

No problem! We have a majority of our available dimensions up on www.phenomenex.com, but if you can't find what you need right away, our super helpful Technical Specialists can guide you to the solution via our online chat portal www.phenomenex.com/Chat.

Australia

t: +61 (0)2-9428-6444
auinfo@phenomenex.com

Austria

t: +43 (0)1-319-1301
anfrage@phenomenex.com

Belgium

t: +32 (0)2 503 4015 (French)
t: +32 (0)2 511 8666 (Dutch)
beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681
info@phenomenex.com

China

t: +86 400-606-8099
cninfo@phenomenex.com

Czech Republic

t: +420 272 017 077
cz-info@phenomenex.com

Denmark

t: +45 4824 8048
nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063
nordicinfo@phenomenex.com

France

t: +33 (0)1 30 09 21 10
franceinfo@phenomenex.com

Germany

t: +49 (0)6021-58830-0
anfrage@phenomenex.com

Hong Kong

t: +852 6012 8162
hkinfo@phenomenex.com

India

t: +91 (0)40-3012 2400
indiainfo@phenomenex.com

Indonesia

t: +62 21 5019 9707
indoinfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405
eireinfo@phenomenex.com

Italy

t: +39 051 6327511
italiainfo@phenomenex.com

Japan

t: +81 (0) 120-149-262
jpinfo@phenomenex.com

Luxembourg

t: +31 (0)30-2418700
nlinfo@phenomenex.com

Mexico

t: 01-800-844-5226
tecnicomx@phenomenex.com

The Netherlands

t: +31 (0)30-2418700
nlinfo@phenomenex.com

New Zealand

t: +64 (0)9-4780951
nzinfo@phenomenex.com

Norway

t: +47 810 02 005
nordicinfo@phenomenex.com

Poland

t: +48 22 104 21 72
pl-info@phenomenex.com

Portugal

t: +351 221 450 488
ptinfo@phenomenex.com

Singapore

t: +65 6559 4364
sginfo@phenomenex.com

Slovakia

t: +420 272 017 077
sk-info@phenomenex.com

Spain

t: +34 91-413-8613
espinfo@phenomenex.com

Sweden

t: +46 (0)8 611 6950
nordicinfo@phenomenex.com

Switzerland

t: +41 (0)61 692 20 20
swissinfo@phenomenex.com

Taiwan

t: +886 (0) 0801-49-1246
twinfo@phenomenex.com

Thailand

t: +66 (0) 2 566 0287
thaiinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367
ukinfo@phenomenex.com

USA

t: +1 (310) 212-0555
www.phenomenex.com/chat

🌐 **All other countries/regions
Corporate Office USA**

t: +1 (310) 212-0555
www.phenomenex.com/chat

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country/region, contact Phenomenex USA, International Department at international@phenomenex.com

**BE-HAPPY™
GUARANTEE**

Your happiness is our mission. Take 45 days to try our products. If you are not happy, we'll make it right.

www.phenomenex.com/behappy

Terms and Conditions

Subject to Phenomenex Standard Terms and Conditions, which may be viewed at www.phenomenex.com/TermsAndConditions.

Trademarks

Kinetex, Luna, and BE-HAPPY are trademarks of Phenomenex. Orbitrap and Q Exactive are trademarks of Thermo Fisher Scientific. SCIEX is a registered trademark and NanoLC is a trademark of AB SCIEX Pte. Ltd.

Disclaimer

Comparative separations may not be representative of all applications. Phenomenex is in no way affiliated with Thermo Fisher Scientific.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures.

© 2022 Phenomenex, Inc. All rights reserved.

