


Experience the Power of Ultra-High Recoveries + Best-in-Class Specificity

Oligonucleotide Analysis

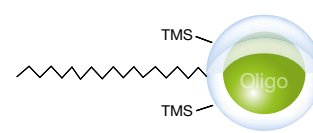

Clarity OTX™ is the industry gold standard for oligonucleotide sample prep

- Ultra-high recoveries for even low nanogram extractions
- Simple 15-minute, 4-step protocol
- Mixed-mode SPE removes salt, sugar protein, and other biological interferences

Proven reliability for analytical characterization of oligonucleotides

- Coreshell morphology improves both mass transfer kinetics and chromatographic efficiency
- BioTi bio-inert titanium hardware mitigates impact of non-specific interactions to improve recovery and peak shape
- Stable from pH 1-12 to accommodate flexible method development demands

DIVERSE SAMPLES

Oligo Types

- DNA
- Aptamers
- RNAi/siRNA
- Thioates
- Lipid-conjugates
- Liposome encapsulated

Sample Types


- Plasma
- Serum
- Urine
- Tears
- Saliva
- Tissue

SAMPLE PREP

The Lysis-Loading buffer was designed with the optimal blend of chaotrope and detergent in order to facilitate the selective binding of oligo therapeutics and metabolites




- Oligo & metabolites
- Salts
- Sugars
- Genomic DNA
- Lipids
- Proteins

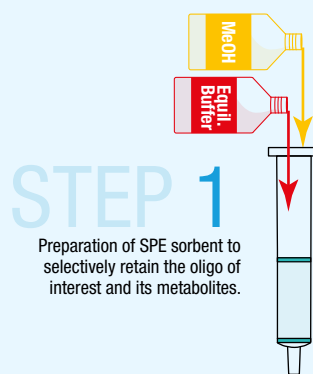


STEP 2

Salts, sugars, large proteins and genomic DNA flow through the cartridge. The oligo of interest, proteins, and lipids bind to the sorbent via a mixed-mode, weak anionic interaction.

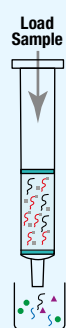


STEP 3

The wash buffer is formulated to strip off lipids and proteins from the sorbent, while not disturbing the oligo therapeutics and its metabolites.

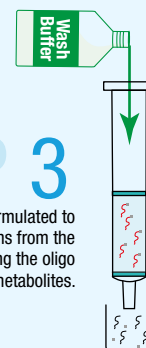


STEP 4

The addition of the Elution buffer releases the target oligo therapeutic and its metabolites. The elution volume can be dried down or lyophilized and reconstituted prior to LC/MS analysis.

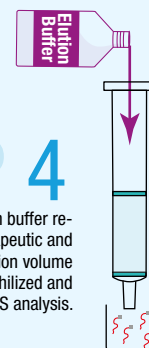


Sensitivity Study in Plasma

Due to high recoveries of parent oligonucleotides and their metabolites, detection down to picomole ranges are possible.

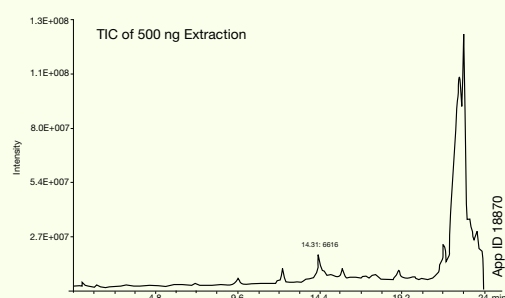


Columns: Clarity 3 µm Oligo-RP C18
Dimensions: 50 x 4.6 mm
Part No.: [00B-4441-EQ](#)
Mobile Phase: A: 50 mM TEAA, pH 7.5 / 5 % Acetonitrile
 B: Methanol
Gradient: A/B (90:10) to A/B (40:60) in 20 min
Flow Rate: 1 mL/min
Detection: UV @ 260 nm
Sample: 25nt DNA Oligonucleotide

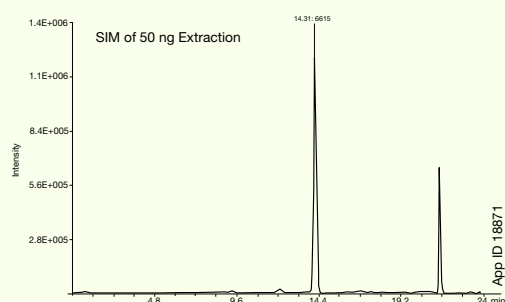

LC-MS Analysis of siRNA using BioTi™ UHPLC Hardware

Bioinert hardware clearly separates sense and anti-sense strands while stainless steel alternatives are susceptible to non-specific interactions that impact reliable quantitation.

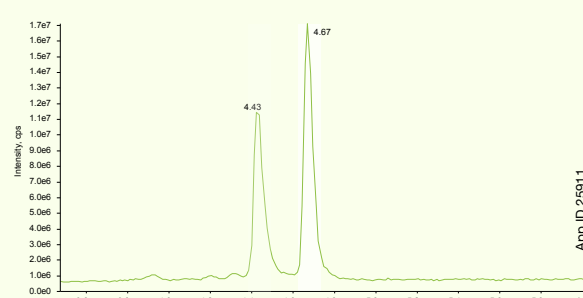

Columns: Biozen 2.6 µm Oligo (BioTi™)
 Clarity 2.6 µm Oligo-XT (stainless steel)
Dimensions: 100 x 2.1 mm
Part No.: [00D-4790-AN](#) (Biozen)
[00D-4746-AN](#) (Clarity)

Mobile Phase: A: 4 mM Triethylamine in Water + 12.5 mM Hexafluoro-2-propanol
 B: 4 mM Triethylamine in Methanol + 12.5 mM Hexafluoro-2-propanol

Gradient:	Time (min)	%B
	5	5
	2	5
	16	90
	16.1	95
	20	95
	20.1	5

Flow Rate: 0.3 mL/min

Injection Volume: 2 µL (12.5 ng)

Temperature: 55°C

Instrument: Shimadzu® LC-20A Prominence®

Detection: TOF-MS

Detector: SCIEX™ TripleTOP® 6600

ANALYSIS

Clarity OTX SPE Ordering Information

Part No.	Description	Unit
KS0-8494	Clarity OTX Starter Kit-Tubes Includes: 100 mg/3 mL cartridges (x50) Lysis-loading buffer (100 mL) Equilibration buffer (250 mL) Wash buffer (350 mL) Elution buffer (100 mL)	ea
KS0-9253	Clarity OTX Starter Kit-96-Well Plate 100 mg/ 96-well plate (x1) Lysis-loading buffer (100 mL) Equilibration buffer (250 mL) Wash buffer (350 mL) Elution buffer (100 mL)	ea
8M-S103-4GA	Clarity OTX Microelution Well Plate	2mg/ well 1/box
8E-S103-CGA	Clarity OTX Well Plate	25 mg/ well 1/box
8E-S103-EGA	Clarity OTX Well Plate	100 mg/ well 1/box
8B-S103-EBJ	Clarity OTX Cartridge	100 mg/3 mL 50/box
8B-S103-HCH	Clarity OTX Cartridge	500 mg/6 mL 30/box
AL0-8579	Clarity OTX Lysis-Loading Buffer V2.0	1 L ea



Biozen Oligo LC Ordering Information

	50 x 2.1	100 x 2.1	150 x 2.1
Biozen 1.7 µm Oligo	00B-4791-AN	00D-4791-AN	00F-4791-AN
Biozen 2.6 µm Oligo	00B-4790-AN	00D-4790-AN	00F-4790-AN

	50 x 4.6	100 x 4.6	150 x 4.6
Biozen 1.7 µm Oligo	—	—	—
Biozen 2.6 µm Oligo	00B-4790-EQ	00D-4790-EQ	00F-4790-EQ

	for 2.1 mm /3pk	for 4.6 mm /3pk	Holder ea
Biozen 1.7 µm Oligo	AJ0-9820	AJ0-9822	KJ0-9000
Biozen 2.6 µm Oligo	AJ0-9820	AJ0-9822	KJ0-9000



Clarity and Biozen are trademarks of Phenomenex. Clarity, OTX, Biozen, and BioTi are trademarks of Phenomenex. Clarity Oligo-XT is patented by Phenomenex. U.S. Patent No. 7,563,367 and 8,658,038 and foreign counterparts. Clarity OTX is patented by Phenomenex. U.S. Patent No. 7,119,145.

FOR RESEARCH USE ONLY. Not for use in clinical diagnostic procedures

© 2023 Phenomenex, Inc. All rights reserved.



POS4980823_W